

# 12962A, 12963A, 12964A room temperature sample holders



# 12962A, 12963A, 12964A sample holders

#### 12962A room temperature sample holder

The 12962A sample holder is designed to allow accurate impedance tests of solid materials to be performed at room temperature. The sample holder consists of two parallel electrodes, one of which is fixed in position and the other which can be moved into contact with the sample by adjustment of a micrometer. The sample holder makes use of guard ring and reference techniques in order to improve accuracy by reducing "fringing" effects at the edge of the sample. The standard electrode provided is 20mm diameter.

#### 12963A electrode kit

Auxiliary electrode kit for the 12962A sample holder allows measurements of different sizes of materials. 10, 30 and 40mm electrodes are supplied in this kit.

### 12964A liquid sample holder

The liquid sample holder also makes use of guard ring techniques and is shaped as a container in order to be able to measure small samples of liquid or powders. The sample holder is easily dissembled for cleaning.





#### **Specification**

electrode diameter :10, 20, 30 and 40mmconnections:2-terminalsample types:solid, liquid and powdersample thickness:0.2mm to 25.4mmsample width and length:≥10mmImpedance range:10hm to 100Tohms

This sample holder is designed for room temperature measurements only, other sample holders and systems are available for high and low temperature measurements in cryostats and furnaces.



Figure 1: Electrode arrangement of he 12962A sample holder

## **Measurement Techniques**

#### Guard ring

The 12962A, 12963A and 12964A sample holders make use of a guard ring on the fixed electrode in order to reduce the effect of stray field lines at the edge of the sample which would otherwise lead to measurement errors. The guard ring ensures that the electric field lines are parallel throughout the part of the sample which contributes to the impedance measurement.

#### **Normalised Measurements**

In order to obtain best accuracy from any sample holder, it is advisable to make use of normalisation measurement techniques. The sample holder allows normalised measurements to be performed with ease. Typically measurements are taken with the sample material in the cell and then repeated without the sample (using an airgap). The air-gap can easily be set up using the micrometer. The empty cell measurements can then be used to normalise the sample results to further reduce errors due to fringing, connections and instrumentation.



Unit B1 Armstrong Mall Southwood Business Park Farnborough GU14 0NR United Kingdom Tel: +44 (0) 1252 556 800 Fax: +44 (0) 1252 556 899 Solartron Analytical's Quality System is approved to BS EN ISO 9001:2008

![](_page_1_Picture_22.jpeg)

...part of **AMETEK**<sup>®</sup> Advanced Measurement Technology

801 South Illinois Avenue Oak Ridge TN 37831 USA Tel: (1) 865-425-1360 Fax: (1) 865-425-2410

solartron.info@ametek.com