

# Application Note H-1



## Subject: A Table of Selected Half-Wave Potentials for Inorganic Substances

### KEY:

All potentials are referenced to the Saturated Calomel Electrode.

\* - Potentials which have an asterisk are differential pulse peak potentials. All other potentials are half-wave potentials.

[ ] – Where a working electrode other than the Dropping Mercury Electrode is used, the electrode material is indicated by brackets.

( ) – Potential values in parentheses indicate that the electrode reaction is an oxidation.

**Acetate (4.5)** – Numbers in parentheses next to a supporting electrolyte are the pH values for that electrolyte.

Abbreviated electrolytes are explained following the table.

ION	SUPPORTING ELECTROLYTES			
<b>Ag (I)</b>	0.1M KNO <sub>3</sub> : 0.10 [Pt]	NH <sub>3</sub> -NH <sub>4</sub> Cl: -0.24		
<b>Al (III)</b>	SVRS-Ac (4.5): -0.46*	0.1M TMAC: -0.75		
<b>As (III)</b>	1M HCl: -0.42*, -0.84*	1M H <sub>2</sub> SO <sub>4</sub> : -0.43, -0.81	1M NaOH: -0.3	H <sub>2</sub> SO <sub>4</sub> -NaCl: -0.20
<b>As (V)</b>	HClO <sub>4</sub> -Pyrogallol: -0.11			
<b>Au (I)</b>	0.1M KOH: -0.1.16			
<b>Au (III)</b>	1M HCl: 0.37 [GCE]			
<b>Ba (II)</b>	0.1M LiCl: -1.92			
<b>Bi (III)</b>	1M HCl: -0.09	Tartrate(4.4): -0.14	NH <sub>4</sub> Cit(3): -0.19	
<b>Br<sup>-</sup></b>	0.1M KNO <sub>3</sub> : (0.12)	KNO <sub>3</sub> -MeOH: (0.15) [Ag]		
<b>BrO<sub>3</sub><sup>-</sup></b>	H <sub>2</sub> SO <sub>4</sub> - KNO <sub>3</sub> : -0.41	0.1M KCl: -1.78		
<b>Cd (II)</b>	NH <sub>4</sub> Cit(3): -0.63*	Acetate(4.5): -0.65	1M HCl: -0.64	NH <sub>4</sub> Tart(9): -0.59*
<b>Ce (III)</b>	2M K <sub>2</sub> CO <sub>3</sub> : (0.16)			
<b>Ce (IV)</b>	2M K <sub>2</sub> CO <sub>3</sub> : -0.16			
<b>Cl<sup>-</sup></b>	0.1M KNO <sub>3</sub> : (0.25)	KNO <sub>3</sub> -MeOH: (0.28) [Ag]		
<b>ClO<sup>-</sup></b>	0.5M K <sub>2</sub> SO <sub>4</sub> (7): 0.08			
<b>ClO<sub>2</sub><sup>-</sup></b>	1M NaOH: -1.0			
<b>CN<sup>-</sup></b>	Borate(9.75): (-0.27)*	0.1M NaOH: (-0.36)		
<b>Co (II)</b>	NH <sub>3</sub> -NH <sub>4</sub> Cl: -1.30	Py-PyHCl: -1.06	5M CaCl <sub>2</sub> : -0.82	NH <sub>4</sub> Cit(9): -1.39
<b>Cr (III)</b>	KSCN(3.2): -0.75*			
<b>Cr (VI)</b>	NH <sub>3</sub> -NH <sub>4</sub> Cl: -0.30*	1M NaOH: -0.84*	NH <sub>4</sub> Tart(9): -0.24	
<b>Cu (I)</b>	NH <sub>3</sub> -NH <sub>4</sub> Cl: (-0.22), -0.50			
<b>Cu (II)</b>	NH <sub>4</sub> Cit(3): -0.06*	Acetate(4.5): -0.07	1M HCl: -0.22	NH <sub>4</sub> Tart(9): -0.36*
<b>Eu (III)</b>	0.1M NH <sub>4</sub> Cl: -0.67			

<b>Fe (II)</b>	Oxalate(4): (-0.23)*	Na <sub>4</sub> P <sub>3</sub> O <sub>7</sub> (9): (-0.37)*		
<b>Fe (III)</b>	Oxalate(4): -0.23*	NaOH-TEA: -1.01*	NH <sub>4</sub> Tart(9): -1.45*	Na <sub>4</sub> P <sub>3</sub> O <sub>7</sub> (9): -0.99
<b>Ga (III)</b>	1M NaSCN(2): -0.83			
<b>Ge (II)</b>	6M HCl: -0.45			
<b>Hg (II)</b>	1M HCl: 0.44 [Au]			
<b>H<sub>2</sub>O<sub>2</sub></b>	PO <sub>4</sub> -Cit(7): (0.18), -1.0			
<b>I<sup>-</sup></b>	0.1M KNO <sub>3</sub> : (-0.03)	KNO <sub>3</sub> -MeOH: (0.0) [Ag]		
<b>In (III)</b>	Acetate(4.5): -0.71	1M HCl: -0.56		
<b>IO<sub>3</sub><sup>-</sup></b>	Phosphate(6.4): -0.79	1M KCl: -1.16		
<b>IO<sub>4</sub><sup>-</sup></b>	K <sub>2</sub> SO <sub>4</sub> - H <sub>2</sub> SO <sub>4</sub> : -0.12			
<b>Ir (IV)</b>	1M HCl: 0.65 [GCE]			
<b>K (I)</b>	0.1M TBAOH: -2.14			
<b>Li (I)</b>	0.1M TBAOH: -2.33			
<b>Mn (II)</b>	NH <sub>3</sub> -NH <sub>4</sub> Cl: -1.66	NH <sub>4</sub> Tart(9): -1.55*		
<b>Mo (VI)</b>	0.3M HCl: -0.26, -0.63	H <sub>3</sub> Cit: 0.04, -0.44		
<b>Na (I)</b>	0.1M TBAOH: -2.12			
<b>Nb (V)</b>	8M HCl: -0.46, -0.70			
<b>Ni (II)</b>	NH <sub>3</sub> -NH <sub>4</sub> Cl: -1.10	NH <sub>4</sub> Tart(9): -0.98*	Py-PyHCl: -0.75	
<b>NH<sub>3</sub>OH</b>	1M NaOH: (-0.43)			
<b>N<sub>3</sub><sup>-</sup></b>	0.1M KNO <sub>3</sub> : (0.25)			
<b>NO<sub>2</sub><sup>-</sup></b>	DPA-SCN(1): -0.54*	2M Cit(2.5): -1.06*	U(VI)-Ac-KCl(2): -0.98*	
<b>NO<sub>3</sub><sup>-</sup></b>	U(VI)-Ac-KCl(2): -0.54*			
<b>O<sub>2</sub></b>	0.1M KNO <sub>3</sub> : -0.05, -0.90			
<b>Pb (II)</b>	NH <sub>4</sub> Cit(3): -0.48*	Acetate(4.5): -0.50	1M HCl: -0.44	NH <sub>4</sub> Tart(9): -0.52*
<b>Pd (II)</b>	NH <sub>3</sub> -NH <sub>4</sub> Cl: -0.75			
<b>Rb (I)</b>	0.1M TBAOH: -2.03			
<b>Rh (III)</b>	NH <sub>3</sub> -NH <sub>4</sub> Cl: -0.93			
<b>Ru (IV)</b>	1M HClO <sub>4</sub> : 0, 0.02, -0.34			
<b>S<sup>2-</sup></b>	0.1M NaOH: (-0.78)*			
<b>Sb (III)</b>	6M HCl: -0.23*	1M HCl: -0.15		
<b>Sb (V)</b>	6M HCl: -0.23*			
<b>Se (IV)</b>	1M HCl: -0.10, -0.40			
<b>Sn (II)</b>	1M HCl: (-0.1), -0.47	NH <sub>4</sub> Cit(3): (-0.21), -0.54	NH <sub>3</sub> Tart(9): -(0.53), -0.77	1M NaOH: (-0.73), -1.22
<b>Sn (IV)</b>	HCl-NH <sub>4</sub> Cl: -0.25, -0.52	1M HCl: -0.1, -0.47		
<b>SO<sub>3</sub><sup>2-</sup></b>	Acetate(5): (-0.62)*			
<b>S<sub>2</sub>O<sub>3</sub><sup>2-</sup></b>	Acetate(5): (-0.21)*			
<b>Te (IV)</b>	NH <sub>4</sub> Tart(9): -0.71	NH <sub>3</sub> -NH <sub>4</sub> Cl: -0.67		
<b>Ti (III)</b>	H <sub>2</sub> Tart: (-0.44)			
<b>Ti (IV)</b>	0.1M HCl: (-0.81)			
<b>Tl (I)</b>	Acetate(4.5): -0.47	1M HCl: -0.48	Ac-EDTA(4.5): -0.50	
<b>U (VI)</b>	0.1M HCl: -0.18, -0.94			
<b>V (V)</b>	H <sub>2</sub> SO <sub>4</sub> -KSCN: -0.52*			
<b>W (VI)</b>	10M HCl: 0, -0.60			
<b>Zn (II)</b>	NH <sub>4</sub> Cit(3): -1.05*	Acetate(4.5): -1.1	NH <sub>3</sub> -NH <sub>4</sub> Cl: -1.35	NH <sub>4</sub> Tart(9): -1.24*

Supporting Electrolytes are listed in approximate order of preference. These supporting electrolytes were selected from commonly available literature sources are not meant to be all-inclusive.

## SUPPORTING ELECTROLYTES

The list below includes only those supporting electrolytes which were abbreviated for the table.

1. Ac-EDTA(4.5): 0.1M Sodium Acetate - 0.1M Acetic Acid – 0.1M Disodium Ethylenediamine tetraacetic acid, Ph 4.5
2. Acetate(4.5): 0.1M Sodium Acetate – 0.1M Acetic Acid, pH 4.5
3. Acetate(5): 0.1M Sodium Acetate + Acetic Acid to pH 5
4. Borate(9.75): 0.1M Boric Acid + NaOH to pH 9.75
5. 2M Cit(2.5): 2M Citric Acid + NaOH to pH 2.5
6. DPA-SCN(1):  $1.3 \times 10^{-4}$  M Diphenylamine (DPA) – 0.01M NaSCN – 0.04M HClO<sub>4</sub>
7. H<sub>3</sub>Cit: Saturated Citric Acid
8. HCl-NH<sub>4</sub>Cl: 1M HCl – 4M NH<sub>4</sub>Cl
9. HClO<sub>4</sub>-Pyrogallol: 2M HClO<sub>4</sub> – 0.5M Pyrogallol
10. H<sub>2</sub>SO<sub>4</sub>-KNO<sub>3</sub>: 0.1M H<sub>2</sub>SO<sub>4</sub> – 0.2M KNO<sub>3</sub>
11. H<sub>2</sub>SO<sub>4</sub>-KSCN: 0.1M H<sub>2</sub>SO<sub>4</sub> – 0.1M KSCN
12. H<sub>2</sub>SO<sub>4</sub>-NaCl: 2M H<sub>2</sub>SO<sub>4</sub> – 2M NaCl
13. H<sub>2</sub>Tart: Saturated Tartaric Acid
14. KNO<sub>3</sub>-MeOH: 0.1M KNO<sub>3</sub> in 50% MeOH
15. KSCN(3.2): 0.2M NaSCN – 0.2M Acetic Acid, pH 3.2
16. K<sub>2</sub>SO<sub>4</sub>- H<sub>2</sub>SO<sub>4</sub>: 0.16M K<sub>2</sub>SO<sub>4</sub> – 1M H<sub>2</sub>SO<sub>4</sub>
17. NaOH-TEA: 0.3 Triethanolamine 0.2M NaOH
18. Na<sub>4</sub>P<sub>3</sub>O<sub>7</sub>(9): 0.2M Sodium Pyrophosphate + H<sub>3</sub>PO<sub>4</sub> to pH 9
19. NH<sub>3</sub>-NH<sub>4</sub>Cl: 1M NH<sub>3</sub> – 0.1M NH<sub>4</sub>Cl
20. NH<sub>4</sub>Cit(3): 0.1M Citric Acid + NH<sub>4</sub>OH to pH 3
21. NH<sub>4</sub>Cit(9): 0.1M Citric Acid + NH<sub>4</sub>OH to pH 9
22. NH<sub>4</sub>Tart(9): 0.1M Tartaric Acid + NH<sub>4</sub>OH to pH 9
23. Oxalate(4): 0.1M Oxalic Acid + NaOH to pH 4
24. Phosphate(6.4): 0.2M Sodium Dihydrogen Phosphate + NaOH to pH 6.4
25. PO<sub>4</sub>-Cit(7): 0.1M Sodium Dihydrogen Phosphate – 0.1M Sodium Citrate adjusted to pH 7
26. Py-PyHCl: 0.1M Pyridine – 0.1M Pyridine-HCl
27. SVRS-Ac(4.5): 0.1M Acetate Buffer, pH 4.7 –  $1.4 \times 10^{-6}$  M Solochrome Violet RS – 12% Ethanol
28. Tartrate(4.4): 0.1M Na<sub>2</sub>Tartrate, pH 4.4
29. 0.1M TBAOH: 0.1M Tetramethylammonium hydroxide
30. 0.1M TMAC: 0.1M Tetramethylammonium chloride
31. U(VI)-Ac-KCl(2): 20 ppm U(VI) – 0.2M KCl – 0.1M Acetic Acid

## REFERENCES

1. Princeton Applied Research Application Briefs and Application Notes.
2. D. T. Sawyer and J. L. Roberts, Jr., "Experimental Electrochemistry for Chemists", John Wiley and Sons, NY, 1974.
3. L. Meites, "Polarographic Techniques", Interscience Publishers, NY, 1955.
4. L. Meites, "Handbook of Analytical Chemistry", McGraw-Hill Book Co., NY, 1963.