

EIS cell analysis techniques State of charge analysis

John Harper and Brian Sayers
Solartron Analytical, Farnborough, UK.

Application Guide: AGML02

Introduction

The use of impedance measurements to determine the state of charge of a battery is well known. Tests are often run to determine the battery impedance in its fully charged, partially charged and fully discharged state. Impedance measurements not only give information about the state of charge of a cell but also gives a lot of information about the cell's state of health. Failure modes can be identified well before actual failure takes place by examining the trend of impedance after various numbers of charge / discharge cycles.

The following experiment can be customized by the user in many ways, for example:

- If impedance is to be checked as a function of the number of charge / discharge cycles run on a particular cell, this can easily be done by adding a repeat loop to the sequence below.
- Impedance tests can be run on fuel cells at various load currents. In this case only the discharge part of the sequence is used and a DC current is entered in the DC settings of the impedance step. Impedance can be checked at different DC current loading levels by adding more impedance steps into the sequence, each with different DC current load setting.
- Alternatively, additional discharge and impedance steps may be added to break up the discharge process into segments, allowing a detailed impedance vs state of discharge map to be built for a particular cell. This gives a great deal of information about how various cells compare during the discharge process.

The ModuLab system is totally flexible and can be used to run virtually any type of impedance test.

Key system capabilities used in this demonstration

- Integration of DC charge / discharge cycling with impedance techniques

Equipment required for this demonstration

- ModuLab electrochemical test system with FRA option fitted
- Sealed lead acid battery - e.g. 6 V or 12 V (for example 2.5 Ah)

Connections

- Connect the ModuLab potentiostat to the battery following the connection diagram shown in the following experiment.

Experiment setup

Select "AGML02 State of Charge" in the "ModuLab Application Guide" project

Step #	Purpose
Step 1	Charge the battery until it reaches a particular voltage
Step 2	Open circuit to allow the cells to settle after charge
Step 3	Run impedance sweep from 100 kHz to 1 Hz measuring main and auxiliary channels
Step 4	Discharge the battery until it reaches a particular voltage
Step 5	Open circuit to allow the cells to settle after charge
Step 6	Run impedance sweep from 100 kHz to 1 Hz measuring main and auxiliary channels

Additional test possibilities:

- Additional AC levels can be tried
- Battery can be discharged to different levels before running the impedance tests


Notes on setup

Using the experimental setups described in the tables below, it is relatively simple to demonstrate the use of impedance for analyzing the state of charge of primary (single use) or secondary (rechargeable cells). In the case of primary cells or fuel cells, of course, the charge step should be removed from the sequence.

The cell is only partially charged and discharged to keep the time duration of the demonstration to a reasonably short period. In practice, charge and discharge times would be much longer and at higher rate, which would more appreciably change the state of charge of the cell.

Data presentation and analysis

The impedance of the cell when charged and discharged is shown above in Bode and complex plane format (graph format can be selected by right clicking on the graph and changing the graph type in the setup menu). The cell is clearly higher impedance when discharged than when it is charged (there is a 10 % difference in cell impedance even on this very short duration partial discharge test) and it can therefore be seen that impedance can readily be used as an indicator of state of charge of the cell. Additional cell "state of health" information is also available from detailed analysis of the shape of the curve.

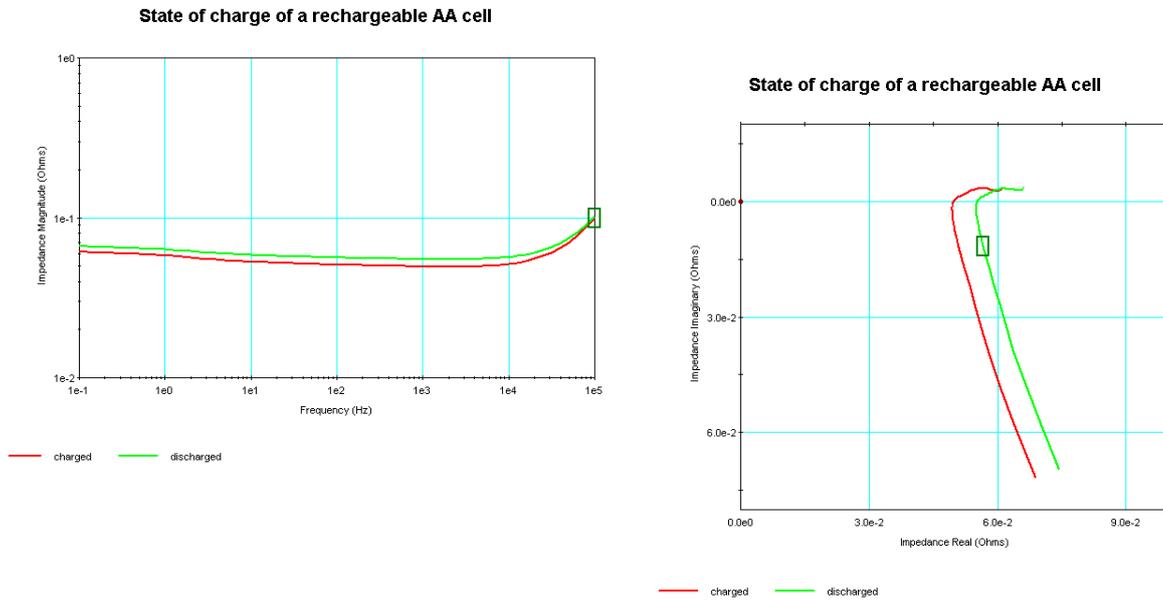

Conclusions

Impedance measurements may be used to determine the state of charge and state of health of a cell (whether it is a single-use or rechargeable cell) and this provides essential information to the engineer or scientist regarding mechanisms that are critical toward the development and realization of emerging technologies.



Solartron Analytical's Quality System is
approved to
BS EN ISO 9001:1994



FM01709

...part of **AMETEK**® Advanced Measurement Technology

UNIT B1 ARMSTRONG MALL
SOUTHWOOD BUSINESS PARK
FARNBOROUGH, GU14 0NR
UNITED KINGDOM
Phone: +44 (0) 1252 556 800
Fax: +44 (0) 1252 556 899

801 SOUTH ILLINOIS AVENUE
OAK RIDGE
TN 37831-2011
USA
Phone: +1 865 425 1360
Fax: +1 865 425 1334

Visit our website for a complete list of our global offices and authorized agents

solartron.info@ametek.com

www.solartronanalytical.com