

How To Guide



Installing the G0159 Reference Electrode Jacket in the 303A



To install the G0159 Reference Electrode Jacket into the 303A cell block, it is best to first wet the O-ring then, while applying gentle upward pressure, use a small flat blade screw driver or other instrument to work the O-ring into the opening so as to wedge the assembly into the opening.



When installed properly, the O-ring should be fully inserted into the opening as shown above. To remove the G0159 assembly, grasp the reference jacket and apply downward pressure while working it back/forth. This action should be sufficient to remove it for maintenance.